Rev. 2026/03

Instant FLYFood

Instant Powdered Medium for *Drosophila melanogaster* – Biosigma

Description:

Instant FLYFood is a powdered culture medium designed for the rapid and standardized preparation of nutritive medium for the rearing of *Drosophila melanogaster*.


Ready to use-mixture packaged in a durable plastic bottle. Simply weight the required amount and add room-temperature or slightly warm water.

The specialized stabilized formulation provides:

- Faster preparation compared to traditional methods, without the need of boiling
- Reduced operational costs through lower energy consumption and simplified equipment requirements
- Compatibility with thermosensitive compounds, which can be directly added to the hydration water
- Natural control of microbial growth without additional supplements
- High reproducibility across preparations and production batches

Instructions for Use:

1. Weigh 0.192 g of powder per 1 mL of water directly into the final container (vial or bottle).
2. Place the container on a vortex mixer and start agitation.
3. Add water while **maintaining agitation** to ensure homogeneous dispersion of the powder.
4. Allow to stand until full water absorption and uniform gel formation are achieved (about 45-60 minutes).
5. Use or store preparations not intended for immediate use in a refrigerator.

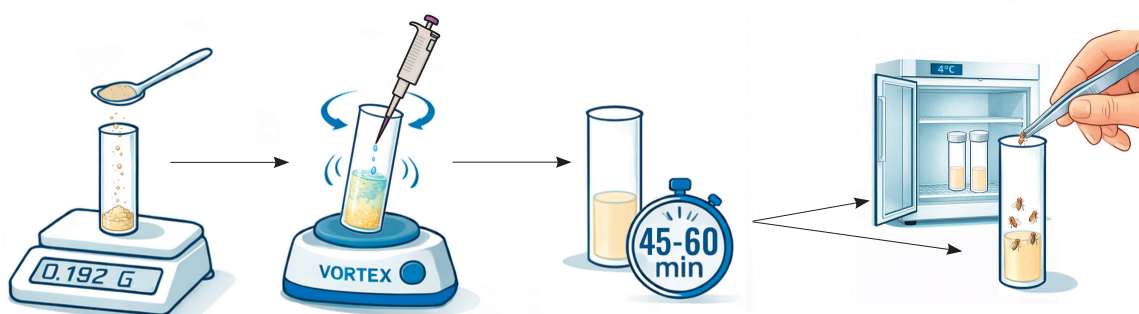Technical Note:

To avoid dispensing difficulties due to increased viscosity during gel formation, it is recommended to:

- Weigh the powder directly into the final container otherwise
- Work with small volumes under continuous agitation and dispense the preparation before complete gelification.

Storage and Stability:

- **Storage conditions:** Store in a dry place, protected from light and heat sources.
- **Recommended temperature:** 15-25 °C.
- **Shelf-life:** 24 months from the date of manufacture, if stored according to the indicated conditions.

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Biosigma S.p.A Dominique Dutscher Group