Biosigma

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Rev 2026/03

Instant FLYFood - Ready-to-Use Vials

Pre-dosed Instant Medium fo *Drosophila melanogaster* - Biosigma

Description:

Instant FlyFood - Vials is the first instant powdered medium, pre-dosed directly into vials, designed for the rapid and standardized preparation of nutritive medium for *Drosophila melanogaster* rearing. The mixture, pre-dosed and packaged directly in the final rearing vial, only requires the addition of room-temperature or slightly warm water to form a stable and homogeneous gel.

The specially stabilized formulation ensure:

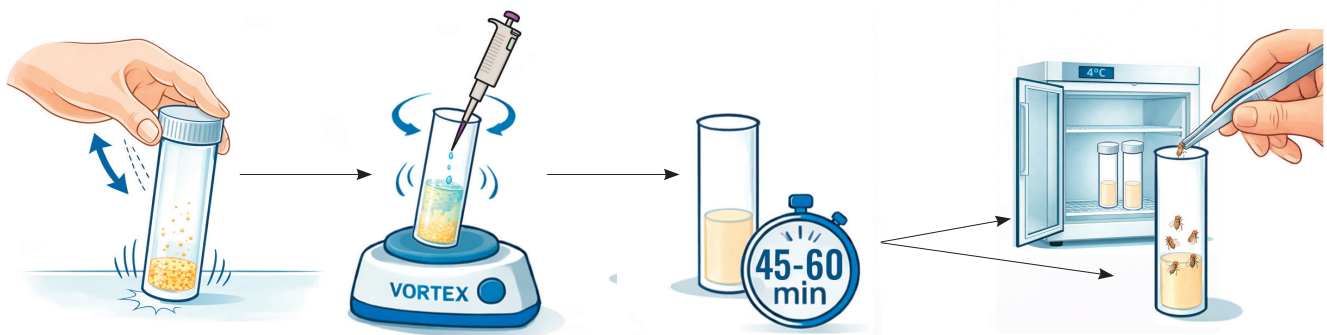
- Immediate preparation compared to traditional methods, with no weighing, transferring, or boiling required
- Reduced preparation time and operational costs
- Compatibility with thermosensitive compounds, which can be directly added to the hydration water
- Natural control of microbial growth, without the need for additional supplements
- High reproducibility and standardization between vials and between batches

The product is available in the following formats:

- Ø 25 mm - Code BFF025I
- Ø 28 mm - Code BFF028I

Instructions for Use:

1. Gently tap the vial to allow the powder to settle completely at the bottom, then remove the cap.
2. Place the vial on a vortex mixer and start agitation.
3. While **maintaining agitation**, add:
 - 5 mL of water for Ø 25 mm vials
 - 8 mL of water for Ø 28 mm vials
4. Vortex vigorously until the powder is completely and uniformly hydrated.
5. Allow to stand until full water absorption and uniform gel formation are achieved (about 45-60 minutes).
6. The product is ready for *Drosophila* rearing: apply the suitable cap (not supplied) for immediate use or store the vials in a refrigerator until use.

Storage and Stability:

- **Unhydrated product:** Store in a dry place at 15–25 °C, protected from light and heat sources.
- **Hydrated product:** Store at 4 °C until use.
- **Shelf-life:** 24 months for the unhydrated product when stored under the indicated conditions.

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Biosigma S.p.A Dominique Dutscher Group